Bottle Adapter EBI – FLS

1 Description

The bottle adapter is easy to install and suitable for bottles with a neck diameter of approximately 27 mm .

It is made to attach a logger of the EBI 12-T 23x, EBI 12-T43x or EBI 12-T46x at a bottle.

2 Scope of delivery

Check contents of the packing for completeness and integrity.


If you discover any damage or have any reason for complaint, please contact your distributor or:

Xylem Analytics Germany GmbH
Am Achalaich 11
82362 Weilheim
Germany
E-Mail: ebro@xylem.com
Internet: www.ebro.com

Parts list:

EBI-FLS, Silikon, up to 130 °C


- Accessories:
 - 1 x Nut M 10
 - 10 x O-Ring
 - 1 x Grommet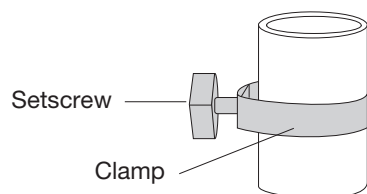

3 Assembling

3.1 Put on the adapter


Put on the adapter on to the bottle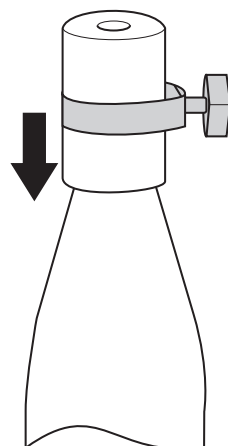

Turn the setscrew to the left, if the diameter of the clip is too small, in order to increase it. Push the adapter as far, as it rests shortly at the bottle neck.

Turn the set screw to the right, in order to fix the adapter in this position.

Tighten the set screw always firmly.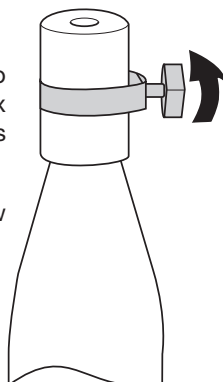

3.2 Preparation of the logger


Screw enclosed M 10 - nut on the thread of the logger.

The straight side of the nut thereby points to the logger, the chamfered side upward.

The position of the nut determines the submergence of the sensor.


Move the O-ring, until it rests shortly to the nut.

Finally attach the grommet.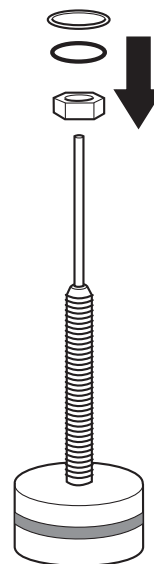

3.3 Assemble the measuring system

Attach the completed logger into the bottle adapter.

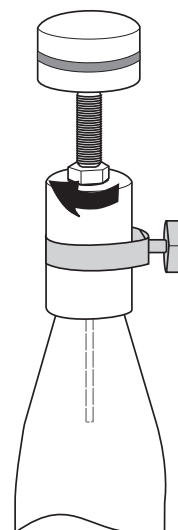Turn the logger to the right into the bottle adapter, until desired submergence is reached.

If necessary turn at the nut, in order to position the sensor exactly.

Screw the nut with a wrench (WS 17) firmly. Now the system is ready to measure.

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